

***ClearBand* WellTracker, 100x**

1 ml, 5x1 ml

Cat No: WT1, WT5

Shipping : Ship at ambient temperature.

Storage : Store at room temperature.

General Information

ClearBand WellTracker is specifically formulated to easily see the wells of the SDS-PAGE gels when loading samples. It is supplied as 100x concentrate and added in stacking gel to better visualize the wells of the gel. The special dye within *ClearBand* WellTracker migrates with bromophenol blue and does not interfere with the samples when they run in the gel. It is totally inert and does not affect the transfer efficiency of proteins from gel to the membrane.

Protocol

1. Add 30 μ l *ClearBand* WellTracker to 3 ml stacking gel just before adding TEMED and APS. Mix well and pour the gel as you routinely do.
2. Wait for solidification of the gel, load your samples and run your SDS-PAGE.