

NutriCulture 100x CellSAFE

50 ml, 100 ml

Cat No: CS50, CS100

Shipping : Ship at ambient temperature.

Storage : Store at room temperature.

General Information

Contamination of cell cultures is a frequent and important problem faced in cell and tissue culture laboratories. It results in delays in experimental schedules, and thus causes waste of precious time, money and efforts. 100x CellSAFE helps keeping your incubators clear of undesired contamination with its strong antimicrobial and fungicidal features against a wide variety of well-known lab contaminants, without effecting the morphology and proliferative potential of routinely used cells in cell and tissue culture laboratories.

Recommended use

1. Cleanse the incubator water pan and fill it with autoclaved distilled water.
2. Add 10 mL of 100x CellSAFE per 1 L of water in the water pan.
3. Replace water within the pan every 2 to 4 weeks and add 100x CellSAFE each time.

Precaution: Note that some precipitate may be observed in the water pan, which can be removed easily with gentle rinsing.