

***ClearBand* Laemmli Sample Buffer (10x)**

2x1 ml

Cat No: LSB-10x

Shipping : Ship at ambient temperature.

Storage : Store at room temperature.

General Information

ClearBand Laemmli Sample Buffer 10x is a premixed buffer for protein sample preparation prior to loading on SDS-PAGE. It is supplied as 10x concentrate. 20 µl B-mercaptoethanol should be added per 80µl *ClearBand* Laemmli Sample Buffer 10x prior to use.

Protocol

1. Add 20µl B-mercaptoethanol per 80µl *ClearBand* Laemmli Sample Buffer 10x prior to use.
2. Mix your protein sample and *ClearBand* Laemmli Sample Buffer 10x in 9:1 ratio.
3. Heat sample - Laemmli buffer mix for 5 minutes in boiling water.
4. Load boiled samples into the wells of SDS-PAGE.